


### BAKERBOND™ PAH-16 PLUS HPLC Column

- Specially designed for the separation of the sixteen PAH's as defined by the U.S. Environmental Protection Agency (EPA)
- Each sorbent batch is tested with the selectivity test mixture SRM 869 to guarantee column-to-column and batch-to-batch reproducibility
- Each column is individually tested with SRM 1647d
- the certificate of analysis and testchromatogram are provided with each column
- the use of narrow column diameter (3 mm) results in better sensitivity

The separation of PAHs is greatly influenced by the type of synthesis that is used for the preparation of C<sub>18</sub>. In reversed phase chromatography two types of silane modification can be distinguished that resulted in either monomeric or polymeric bond linkages. The polymeric phases provided significant enhanced selectivity characteristics, particularly for isomeric PAHs. BAKERBOND™ PAH-16 Plus column contains a wide pore C<sub>18</sub> bonded phase synthesised under polymer-forming conditions. All wide pore C<sub>18</sub> batches are subjected (internal quality control) to a characterization test ( $\alpha_{\text{TBN/BaP}}$ -test) developed by Sander and Wise. The test is intended primarily for predicting the potential selectivity for complex PAH mixtures and to discriminate between monomeric and polymeric type of C<sub>18</sub> stationary phases.

The  $\alpha_{\text{TBN/BaP}}$ -test is used by J.T. Baker to verify column reproducibility and assist in the adjustment of the bonding process. Complete separation of the 16 PAHs, identified by the U.S. EPA as priority pollutants or even more, can only be achieved using bonded phases with  $\alpha_{\text{TBN/BaP}}$  values in a well defined range.

Peak No.	Retention Time (min)
1	4.00
2	4.30
3	4.70
4	5.00
5	5.30
6	5.60
7	5.90
8	6.20
9	6.50
10	6.80
11	7.10
12	7.40
13	7.70
14	8.00
15	8.30
16	8.60

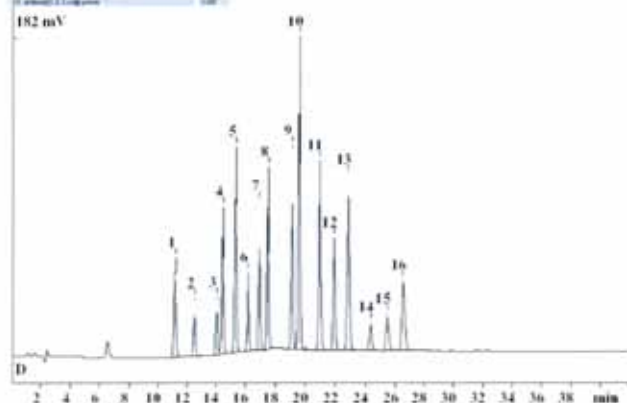

BAKERBOND™ PAH 16 Plus, 5µm,  
250 x 3.0 mm + guard  
**Part No. 7504-00**

Mobile Phase: A=Water  
B=Acetonitrile

Gradient: 50% B for 5 min  
50 – 100% Bin 15 min  
100% B for 25 min

Flow Rate: 0.5 mL/min  
Temperature: 25 °C  
Detection: UV at 254nm

Ref:  
W. Hesselink, R.H.N.A. Schiffer, P. Kootstra, *J. Chromatogr. A*,  
697 (1995) 165 - 174

tyco / Healthcare / Mallinckrodt Baker

www.jtbaker.com

**Mallinckrodt Baker**  
Holland Headquarters  
Teugseweg 20  
PO Box 1  
7400 AA Deventer  
The Netherlands  
t 0031 570 687500  
f 0031 570 687574  
e jtbaker.nl@  
www.tycohealthcare.com

**Mallinckrodt Baker**  
UK  
107112 Leaderhall Street  
London  
EC3A 4AH  
United Kingdom  
t 0044 1908 506000  
f 0044 1908 503290  
e jtbaker.uk@  
www.tycohealthcare.com

**Mallinckrodt Baker**  
Italy  
Via P. Porciacappi 112  
20138 Milano  
Italy  
t 0039 02 580891  
f 0039 02 58013204  
e jtbaker.it@  
www.tycohealthcare.com

**Mallinckrodt Baker**  
Germany  
Im Leuschnerpark 4  
D-64347 Griesheim  
Germany  
t 0049 6155 867939  
f 0049 6155 967940  
e jtbaker.de@  
www.tycohealthcare.com

**Mallinckrodt Baker**  
France  
33 Avenue Philippe Auguste  
75011 Paris  
France  
t 0033 144 64 99 7107  
f 0033 144 64 99 73  
e jtbaker.fr@  
www.tycohealthcare.com



A CENTURY  
OF EMPOWERING  
DISCOVERY™

### PRODUCT LIST

BAKERBOND™ PAH 16 plus	micron	column size (mm)	A	Product number
BAKERBOND™ PAH 16 plus including guard column, manufit connection system and certificate of analysis	5	3 x 250	120	7504-00
Replacement guard cartridges for BAKERBOND™ PAH 16 plus (3)	5	3 x 20	120	7505-00
Replacement column for BAKERBOND™ PAH 16 plus including certificate of analysis	5	3 x 250	120	7506-00



BAKERBOND™ PAH 16/ 251004/SepTec

**tyco** | Healthcare | **Mallinckrodt Baker**

[www.jtbaker.com](http://www.jtbaker.com)

**Mallinckrodt Baker**  
Holland Headquarters  
Teugeweg 20  
PO Box 1  
7400 AA Deventer  
The Netherlands  
t 0031 570 687500  
f 0031 570 687574  
e jtbaker.nl@  
ema@tycohealthcare.com

**Mallinckrodt Baker**  
UK  
107112 Leadenhall Street  
London  
EC3A 4AH  
United Kingdom  
t 0044 1908 506000  
f 0044 1908 503290  
e jtbaker.uk@  
ema@tycohealthcare.com

**Mallinckrodt Baker**  
Italy  
Via P. Portogallo 110  
20138 Milano  
Italy  
t 0039 02 580981  
f 0039 02 58019264  
e jtbaker.it@  
ema@tycohealthcare.com

**Mallinckrodt Baker**  
Germany  
Im Leuschnerpark 4  
D-64347 Gräfenheim  
Germany  
t 0049 6155 667939  
f 0049 6155 667940  
e jtbaker.de@  
ema@tycohealthcare.com

**Mallinckrodt Baker**  
France  
33 Avenue Philippe Auguste  
75011 Paris  
France  
t 0033 144 04 39 7112  
f 0033 144 04 39 73  
e jtbaker.fr@  
ema@tycohealthcare.com



A CENTURY  
OF EMPOWERING  
DISCOVERY™